

INSTRUCTIONS

Chemical Spill Kit (45lt)

1. Unpack chemical spill kit and lay out contents away from direct contact with spill. Dress in PPE supplied.
2. When dealing with a spill, first contain the spill with the 1.2m socks as provided.
3. Place chemical absorbent pads and / or scatter on the spill. Allow to absorb the spill.
4. Once you have completed the clean-up, remove the soiled absorbent pads and / or scatter first before lifting the containment socks, ensuring that the spill is completely absorbed.
5. Dispose of the soiled absorbent products in the disposal bags and seal with ties as provided. Clearly mark the disposal bag identifying the spilled product for disposal.
6. **REMEMBER TO REPLACE USED STOCK FROM KIT**
Herewith the list of contents for your convenience.

Qty	Description
1x	Spill Kit Carry Bag – Clear PVC
1x	30lt Chemical Scatter
5x	200g Chemical Absorbent Pads
2x	1.2m Chemical Absorbent Socks
1x	Pair PVC Gloves
1x	Dust Mask
1x	Anti-mist Goggles
1x	Sparkproof Dustpan and Brush
2x	Heavy Duty Disposal Bags with Ties